



Laboratory Information Management System





IntelliLIMS is a Laboratory Information Management System (LIMS). The software was developed in accordance with FDA regulations and is also suitable for use in the following industries:


- Pharmaceutical • Medical Device • Food and Beverage
- Chemical Manufacturing • Environmental Laboratories • Nutraceutical
- Cosmetics • Automotive • Petrochemical • Mining
- Any Business Requiring Management of Samples and Testing

IntelliLIMS is positioned as a easy to use application, that is flexible and scalable to your business needs.

H&A Scientific, Inc. offers a turn key solution with each IntelliLIMS system. This includes on-site installation, training, support by H&A Scientific, Inc. and SOP templates for use and operation

IntelliLIMS provides all of the following features:



- Sample login
- Sample tracking, including Barcode Support
- Scheduling
- Chain of custody
- Result entry with audit trails
- Multi-level QA/QC/Specification checking
- Result Reporting
- Interface with the PC/Chrom Chromatography Data System
- E-mail notification of samples and more...

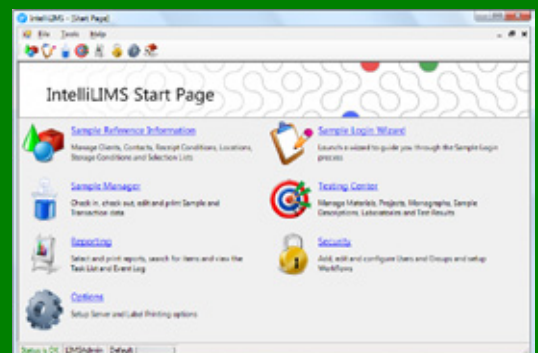

The Intelligent Laboratory Information Management System to Manage Your Laboratory


Additional Features

- IntelliLIMS imports test results from the PC/Chrom Chromatography Data System (CDS).
- The Chromatogram can be viewed directly with the IntelliLIMS Interface.

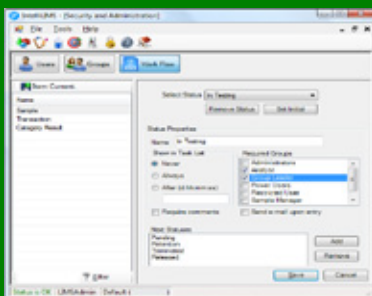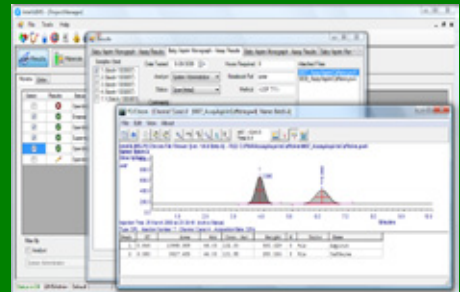

- The workflow is dynamic and configurable.
- E-mail notification can be sent on status changes.
- IntelliLIMS has granular security.
- Windows Authentication is supported.

- Full audit trail including check sums on all records.
- IntelliLIMS includes functionality to meet the Electronic Records and Electronic Signatures rules of 21 CFR Part 11.

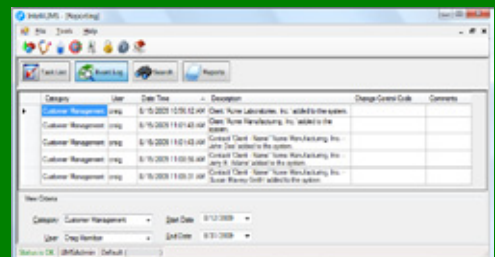


Category	User	Date/Time	Description	Change Control Code	Comments
Customer Management	intg	6/15/2021 10:50:12 AM	Check Name Laboratories, Inc. added to the system.		
Customer Management	intg	6/15/2021 1:01:43 AM	Check Name Pharmaceuticals, Inc. added to the system.		
Customer Management	intg	6/15/2021 1:03:14 AM	Check Client Name Name Manufacturing, Inc. added to the system.		
Customer Management	intg	6/15/2021 1:03:56 AM	Check Client Name Name Manufacturing, Inc. added to the system.		
Customer Management	intg	6/15/2021 1:03:57 AM	Check Client Name Name Manufacturing, Inc. added to the system.		

Validation and Support


- We supply Validation Certificates, Manuals, and Installation Qualification (IQ) documents with all options.
- We also offer additional consulting services in the form of User Acceptance Validation documents and software testing.

Application Flow


Clients and Contacts can be added to IntelliLIMS. This makes it easy to create Client Scoped Projects.

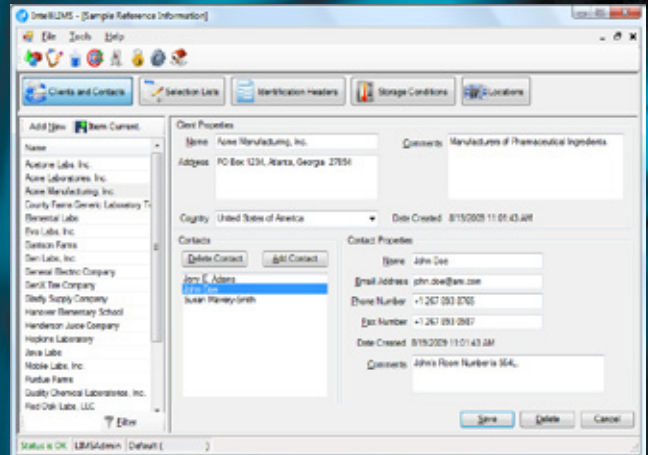

- The Sample Login Wizard allows you to login one or more samples.
- Sample(s) can be assigned a unique identifier.
- Sample Descriptions can be added with associated MSDS link.
- Each Sample can have three ID fields with customizable headings.

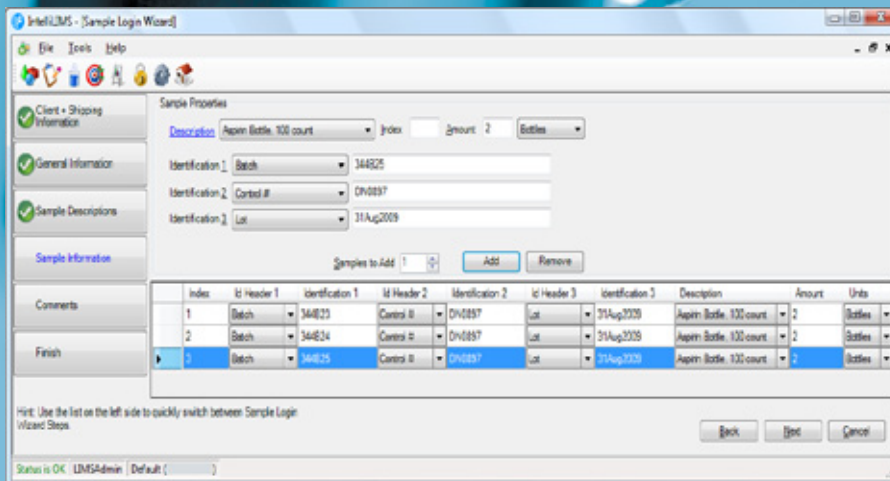


- Barcode labels for samples and transactions can be printed.
- E-mail notifications can be sent.
- Sample Reports can be generated.

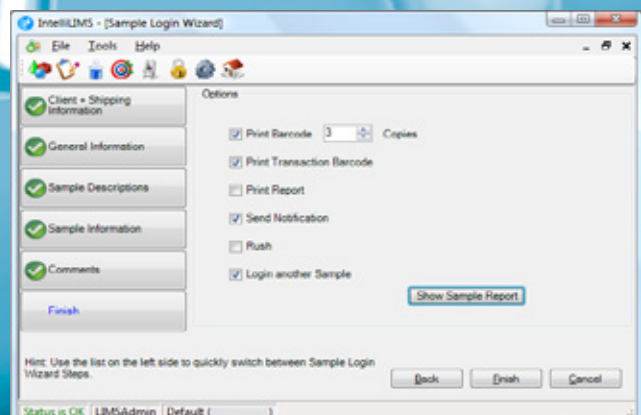


Sample Manager

The Sample Manager maintains control of the samples through the entire workflow process.

- Sample can be easily retrieved by scanning barcodes.
- Checking-in and checking-out samples is supported.
- Each sample has a Chain of Custody
- Sample Reports can be generated.

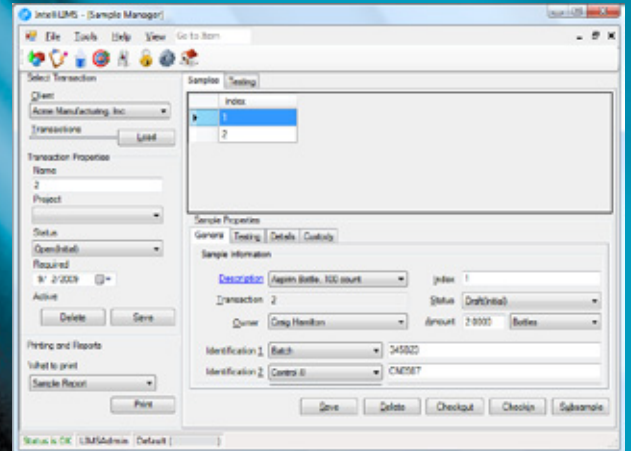

Testing Center

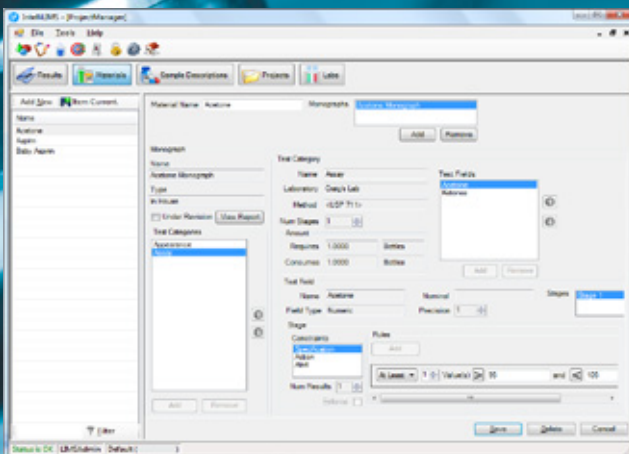

The Testing Center is the foundation for Sample Testing

- Select from a preloaded monograph or create your own custom monograph.
- Both short and long term projects are supported.
- Alert, Action, and Specification limits can be assigned to a test.
- A final Results Summary Report can be generated.

H&A Scientific, Inc. develops, manufactures, and markets quality computing systems for scientific data acquisitions, processing, and management.

H&A Scientific, Inc. is continually maintaining and improving its established Quality Assurance Program which is designed to maintain compliance with requirements of the FDA's Good Manufacturing Practices (GMPs).



For more information, please visit our web site www.hascientific.com or contact:

Daniel Robinson
Manager of Business Development
PO Box 8133, Greenville, NC 27835
Ph: +1.252.752.4315 ext. 204
Fax: +1.252.752.9917
E-mail: dan@hascientific.com

European Support
H&A Scientific, Inc.
Mainzer Landstrasse 176
60327 Frankfurt am Main / Germany
Ph: +49(0)69.97358.261
Fax: +49(0)69.97358.101
E-mail: hainfo@hascientific.com

